



KAPA HiFi™ HotStart ReadyMix (2X)

1. Product Description

KAPA HiFi HotStart ReadyMix (2X) is designed for routine, high-fidelity PCR of a wide range of targets and fragment sizes. It offers error rates 100X lower than that of wild-type *Taq*, as well as higher success rates, yields and consistency – for both AT- and GC-rich amplicons – than those achievable with wild-type B-family (proofreading) DNA polymerases. In addition, KAPA HiFi HotStart ReadyMix reaction times are significantly shorter than those required for wild-type B-family DNA polymerases or hot start formulations thereof.

KAPA HiFi HotStart ReadyMix (2X) is a ready-to-use cocktail containing all components for PCR, except primers and template. The 2X ReadyMix contains KAPA HiFi HotStart DNA Polymerase in a proprietary reaction buffer, dNTPs (0.3 mM of each dNTP at 1X), $MgCl_2$ (2.5 mM at 1X) and stabilizers.

KAPA HiFi HotStart DNA Polymerase is an antibody-based hot start formulation of KAPA HiFi DNA Polymerase, a novel B-family DNA polymerase that exhibits industry-leading performance when compared with other high-fidelity (B-family) DNA polymerases and polymerase blends. KAPA HiFi DNA Polymerase was engineered to have increased affinity for DNA, without the need for accessory protein domains. The intrinsic high processivity of the enzyme results in significant improvements in yield, sensitivity, speed, target length and the ability to amplify difficult amplicons. In the HotStart formulation, the enzyme is combined with a proprietary antibody that inactivates the enzyme until the first denaturation step. This eliminates spurious amplification products resulting from non-specific priming events during reaction setup and initiation, and increases overall reaction efficiency.

KAPA HiFi HotStart DNA Polymerase has 5'-3' polymerase and 3'-5' exonuclease (proofreading) activities, but no 5'-3' exonuclease activity. The strong 3'-5' exonuclease activity results in superior accuracy during DNA amplification. KAPA HiFi HotStart DNA Polymerase boasts the lowest published error rate of all B-family DNA polymerases (1 error per 2.8×10^7 nucleotides incorporated). This fidelity is approximately 100X higher than that of wild-type *Taq* and up to 10X higher than that of other B-family DNA polymerases and polymerase blends. DNA fragments generated with KAPA HiFi HotStart ReadyMix may be used for routine downstream analyses or applications, including restriction enzyme digestion and sequencing. PCR products generated with KAPA HiFi HotStart ReadyMix are blunt-ended, but may be 3'-dA-tailed for cloning into TA cloning vectors.

2. Applications

KAPA HiFi HotStart ReadyMix (2X) is ideally suited for routine, high-fidelity PCR including:

- Amplification of long DNA fragments for conventional sequencing (direct sequencing or sequencing of cloned PCR products).
- Amplification of short AT- and GC-rich DNA fragments for next-generation sequencing (e.g. target enrichment for resequencing).
- Amplification of DNA fragments up to 15 kb from genomic DNA or up to 18 kb from less complex targets (e.g. plasmid or lambda DNA), for cloning and protein expression or genomic characterization.
- Introduction of single or multiple point mutations using site-directed mutagenesis.

For more information on these and other high-fidelity PCR applications, please refer to the KAPA HiFi Application Notes on **Site-Directed Mutagenesis**, **Routine High-Fidelity PCR** and **High-Fidelity GC-rich PCR**.

Kit codes and components

KK2600 Sample kit	KAPA HiFi HotStart ReadyMix (2X) 1 x 0.25 ml (20 x 25 µl rxns)
KK2601 100 reactions	KAPA HiFi HotStart ReadyMix (2X) 1 x 1.25 ml (100 x 25 µl rxns)
KK2602 500 reactions	KAPA HiFi HotStart ReadyMix (2X) 1 x 6.25 ml (500 x 25 µl rxns)

Storage, handling and specifications

Store all components at -20 °C for long-term use. Please refer to Section 6 for full details.

Quick Notes

- KAPA HiFi HotStart ReadyMix contains a novel HotStart DNA Polymerase, engineered for fast and versatile high-fidelity PCR.
- Amplify fragments up to 15 kb from genomic DNA or 18 kb from less complex targets.
- Denature at 98 °C for 20 sec in each cycle.
- Optimal annealing temperatures are higher than in other PCR buffers, and are typically in the range of 60 – 72 °C.
- Use 15 sec extension time per cycle for amplicons ≤1 kb and 30 – 60 sec/kb per cycle for longer amplicons or to improve yields.
- Add 5% DMSO for targets with a GC content >70%.
- For optimal fidelity, always use good quality DNA and the lowest number of cycles.



3. Reaction setup and cycling parameters

The following reaction setup is recommended for KAPA HiFi HotStart ReadyMix reactions:

Component	Final concentration	Volume in a 25 µl reaction ¹
PCR grade water	–	Up to 25.0 µl
2X KAPA HiFi HotStart ReadyMix (contains 2.5 mM Mg ²⁺ at 1X)	1X	12.5 µl
Forward primer (10 µM)	0.3 µM	0.75 µl
Reverse primer (10 µM)	0.3 µM	0.75 µl
DMSO (100%) (for amplicons with a GC content >70%)	5%	1.25 µl
Template DNA ²	As needed	10 – 100 ng for genomic DNA 1 – 10 ng for less complex DNA

¹ For smaller reaction volumes, scale all volumes down proportionately. Reaction volumes >25 µl are not recommended.

² Use 10 ng genomic DNA or 1 ng less complex DNA per 25 µl reaction as a starting point.

The recommended KAPA HiFi HotStart ReadyMix cycling protocol is as follows:

Step	Temperature	Time	Number of cycles
Initial denaturation ¹	95 °C	2 – 5 min	1
Denaturation¹	98 °C	20 sec	15 - 35 ⁴
Primer annealing ²	60 – 75 °C	15 sec	
Extension ³	72 °C	15 – 60 sec/kb	
Final extension	72 °C	1 – 5 min	1
Cooling	4 – 10 °C	HOLD	1

¹ The KAPA HiFi HotStart ReadyMix Buffer has a higher salt concentration than conventional PCR buffers. Since buffer composition affects DNA melting, it is important to denature complex templates and targets with a high GC content sufficiently. **The above denaturation parameters must therefore be strictly adhered to.**

² Due to the high salt concentration of the KAPA HiFi HotStart ReadyMix Buffer, the optimal annealing temperature for a specific primer pair is likely to be different than when used in a conventional PCR buffer. An annealing temperature of 65 °C is recommended as a starting point. Two-step cycling protocols, with a combined annealing/extension temperature in the range of 68 – 75 °C and a combined annealing/extension time of 30 sec/kb may also be used.

³ Use 15 sec per cycle for amplicons ≤1 kb and 30 – 60 sec/kb per cycle for long amplicons or to improve yields.

⁴ ≤25 cycles are recommended for most high-fidelity applications. In cases where very low template concentrations or low reaction efficiency results in low yields, 30 or 35 cycles may be performed to produce sufficient product for downstream analysis or cloning.

For advanced troubleshooting or assistance with reaction setup or optimization, consult the KAPA HiFi FAQs and other web-based technical resources on <http://www.kapabiosystems.com> or e-mail support@kapabiosystems.com.



4. Important parameters

4.1 Denaturation

Due to the higher salt concentration of the KAPA HiFi HotStart ReadyMix buffer, it is important to use the correct denaturation parameters for KAPA HiFi HotStart ReadyMix reactions. Although the HotStart enzyme only requires 30 sec for activation, an initial denaturation time of 2 – 5 min at 95 °C is required to ensure that template DNA is fully denatured before the first primer annealing step. Use 5 min for complex, genomic DNA and GC-rich targets and at least 2 min for less complex templates. In each cycle, denaturation must be performed for 20 sec at 98 °C. Initial denaturation may also be performed at 98 °C, but this may lead to template damage and is not recommended for high-fidelity amplification, especially of long amplicons.

4.2 Primers and annealing temperature

Primer design and quality is important for successful DNA amplification using KAPA HiFi HotStart ReadyMix. Primers should be carefully designed to eliminate the possibility for primer-dimer formation and spurious annealing as far as possible and should have a GC content of 40 – 60%. Primers with a GC content >60% may require a higher denaturation temperature or longer denaturation time. Primers to be used together should have similar theoretical melting temperatures (T_m).

The actual T_m of a given primer may be affected by specific reaction conditions, including the composition of the PCR buffer, DNA concentration, presence of denaturing agents (e.g. DMSO) and nucleotide modifications (e.g. biotin or fluorescent dyes). When using KAPA HiFi HotStart ReadyMix for the first time with a specific primer-template combination, it is recommended that the optimal annealing temperature for the assay be determined empirically in an annealing temperature gradient PCR.

Alternatively, start with an annealing temperature of 65 °C as a first approach. To improve **sensitivity (yields)**, **reduce** the annealing temperature in increments of 1 °C (but not lower than 60 °C). To improve **specificity**, **increase** the annealing temperature in increments of 1 °C (up to 75 °C). For long primers and primer sets with an actual optimal annealing temperature in the range of 68 – 75 °C, use a 2-step cycling protocol with a combined annealing/extension time of 30 sec/kb per cycle.

Always dilute and store primers in a buffered solution (e.g. TE or 10 mM Tris-HCl, pH 8.0 – 8.5) instead of PCR grade water.

4.3 Template DNA

Amplification from low complexity templates, such as lambda or plasmid DNA, is usually easy and should require little optimization. Applications based on low target copy numbers (e.g. when amplifying single-copy genes from genomic templates, or when using cDNA as template) are generally more challenging. For plasmid or phage DNA, 1 – 10 ng template per 25 µl reaction is adequate, whereas up to 100 ng complex genomic DNA or cDNA may be required.

The quality of template DNA may have a significant impact on the outcome of KAPA HiFi HotStart ReadyMix reactions. Degraded, damaged or sheared template DNA is particularly problematic when amplifying targets >1 kb. Always dilute and store DNA in a buffered solution (e.g. TE or Tris-HCl, pH 8.0 – 8.5) instead of PCR grade water, and minimize freeze-thaw cycles to limit degradation and maintain quality. High quality template DNA is essential for high-fidelity amplification.

4.4 Amplicon length

KAPA HiFi HotStart ReadyMix is recommended for the amplification of fragments up to 15 kb from complex, genomic DNA and fragments up to 18 kb from less complex templates. For the efficient amplification of fragments ≥10 kb, higher template concentrations and careful optimization of the final Mg^{2+} concentration (see 4.6) may be required.

4.5 Amplification of GC-rich and other problematic amplicons

For the amplification of fragments with a GC content >70% it is recommended that reactions are supplemented with DMSO, to a final concentration of 5%. Any PCR grade DMSO solution may be used. KAPA HiFi HotStart ReadyMix may also be used in combination with 1X KAPA Enhancer 1 (a proprietary PCR additive supplied in KAPA2G Robust and Robust HotStart PCR Kits), or PCR grade betaine (at a final concentration of 1 M), to improve yields and/or specificity in difficult assays.

If the addition of DMSO, KAPA Enhancer 1 or betaine does not yield satisfactory results, template concentration, Mg^{2+} concentration, annealing temperature and/or extension time may have to be carefully optimized.

4.6 Mg^{2+} concentration

KAPA HiFi HotStart ReadyMix contains Mg^{2+} at a 1X concentration of 2.5 mM, which has been determined to be optimal for most applications. Specific applications, particularly those involving the amplification of fragments >10 kb from complex genomic templates, may require careful optimization of the final Mg^{2+} concentration. KAPA HiFi HotStart ReadyMix reactions may be supplemented with any PCR grade $MgCl_2$ solution. Add 0.25 µl of a 25 mM $MgCl_2$ solution to increase the final Mg^{2+} concentration in a 25 µl reaction by 0.25 mM. The optimal Mg^{2+} concentration for a specific assay may be determined empirically in a $MgCl_2$ gradient PCR.

4.7 TA cloning

DNA fragments generated with KAPA HiFi HotStart ReadyMix may be used directly for blunt-end cloning, or cloning using restriction endonucleases. For TA cloning, 3'-dA-overhangs may be added to blunt-ended DNA fragments in a simple reaction (5 min incubation at 72 °C) containing KAPA Taq DNA Polymerase, 1X KAPA Taq Buffer and dATP (0.2 mM). It is very important to purify KAPA HiFi HotStart ReadyMix PCR products prior to dA-tailing. If this is not done, the 3'-5' exonuclease activity of residual KAPA HiFi HotStart DNA Polymerase will degrade the newly generated dA-overhangs and yield a mixture of dA-tailed and blunt-ended fragments, which is likely to reduce the efficiency of TA cloning.



KAPA HiFi™ HotStart ReadyMix (2X)

5. Troubleshooting

Problem	Possible solutions
Non-specific amplification or smearing	<ul style="list-style-type: none">➤ Increase the annealing temperature by 2 – 5 °C.➤ If an annealing time >15 sec per cycle was used, reduce this to 15 sec.➤ Use less template DNA (but no less than 1 ng genomic DNA/25 µl reaction).➤ Decrease the extension time to 15 sec/kb per cycle.➤ Decrease the number of cycles.➤ Prepare new, good quality template DNA and/or fresh primer stocks.
Low yield	<ul style="list-style-type: none">➤ Decrease the annealing temperature by 2 – 5 °C, but not lower than 60 °C.➤ Use more template (but no more than 100 ng genomic DNA/25 µl reaction.)➤ Increase the extension time to 30 – 60 sec/kb per cycle.➤ Increase the number of cycles.➤ Optimize the final Mg²⁺ concentration in the reaction.

6. Storage, handling and specifications

6.1 Shipping, storage and handling

KAPA HiFi HotStart ReadyMix (2X) PCR Kits are shipped on dry ice or ice packs, depending on the country of destination. Upon receipt, store the entire kit at -20 °C in a constant-temperature freezer. When stored under these conditions and handled correctly, all kit components will retain full activity until the expiry date indicated on the kit.

KAPA HiFi HotStart ReadyMix (2X) contains isostabilizers and may not freeze solidly, even when stored at -20 °C. Nevertheless, always ensure that the ReadyMix is fully thawed and has been vortexed before use.

KAPA HiFi HotStart ReadyMix (2X) may be stored at 4 °C for regular, short-term use (up to 1 month). Provided that it has been handled carefully and not contaminated, the ReadyMix is not expected to be compromised if left (unintentionally) at room temperature for short periods of time (up to 3 days). Long-term storage at room temperature or 4 °C is not recommended. Please note that reagents stored above -20 °C are more prone to degradation when contaminated by the user; storage at such temperatures is therefore at the user's own risk.

6.2 Quality control

KAPA HiFi DNA Polymerase and its proprietary HotStart antibody are extensively purified through the use of multiple chromatography steps. The final formulation contains <2% contaminating protein, as determined in an Agilent Protein 230 Assay. Each batch of ReadyMix is subjected to stringent quality control tests, is free of contaminating exo- and endonuclease activities and meets strict requirements with respect to DNA contamination.

6.3 Product use limitations and licenses

KAPA HiFi HotStart ReadyMix (2X) PCR Kits are developed, designed and sold exclusively for research purposes and *in vitro* use. Neither the product, nor any individual component, has been tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals. Please refer to the MSDS, which is available on request.

Certain applications of this product are covered by patents issued to parties other than Kapa Biosystems and applicable in certain countries. Purchase of this product does not include a license to perform any such applications. Users of this product may therefore be required to obtain a patent license depending upon the particular application and country in which the product is used.

Licensed under U.S. Patent nos. 5,338,671 and 5,587,287 and corresponding patents in other countries.

Boston, Massachusetts, United States

600 West Cummings Park, Suite 5350
Woburn, MA 01801 U.S.A.
Tel: +1 781 497 2933 Fax: +1 781 497 2934

Cape Town, South Africa

2nd Floor, Old Warehouse Building, Black River Park,
Fir Road, Observatory 7925, Cape Town, South Africa
Tel: +27 21 448 8200 Fax: +27 21 448 6503